


PERPETULAB WHEELBARROW SCIENCE TABLE

P7604x34W SERIES TABLES (OAK OR MAPLE)
ASSEMBLY INSTRUCTIONS

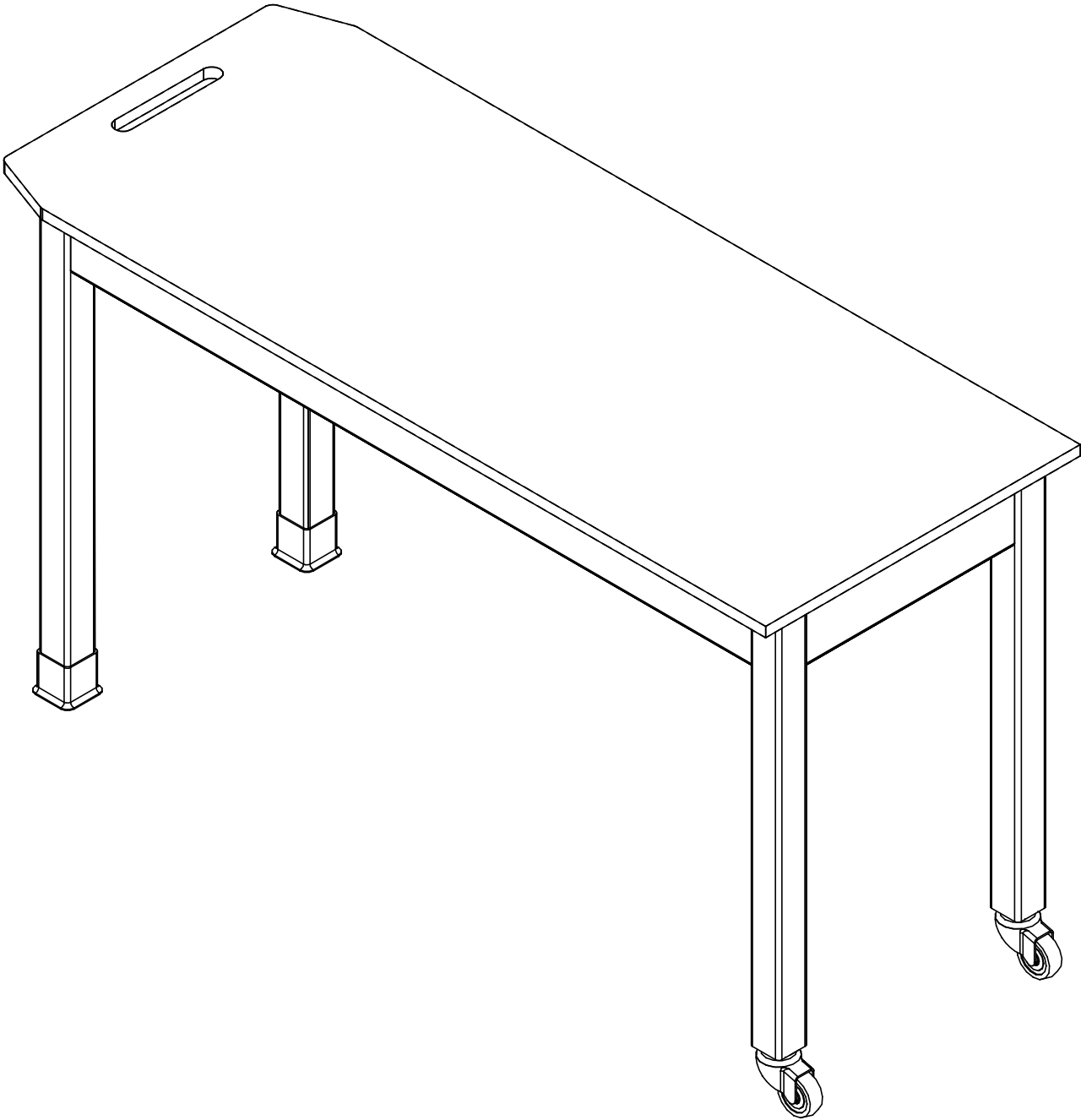

NOTE:

1. MODEL VIEWS MAY NOT REPRESENT EXACT MODEL PURCHASED.

A	CREATED INSTRUCTIONS	09/03/25	EMR	N/A
REV.	DESCRIPTION	DATE	NAME	ECN No.
REVISION LEVEL				

ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE DECIMALS: .X = ± .030 .XX = ± .020 .XXX = ± .010 ANGLES: = ± 1° CABINET SQUARENESS = ± .0625					DWN BY:	EMR	DATE DWN:	09/03/25
					APR BY:		DATE APR:	
					APR BY:		DATE APR:	
CONFIDENTIAL: ALL DESIGN RIGHTS RESERVED THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC. THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.								



TITLE:
PERPETULAB WHEELBARROW SCIENCE TABLE
ASSEMBLY INSTRUCTIONS

DRAWING NUMBER:	SCALE:	SHEET:	REVISION:
P7604x34W SERIES TABLES	NTS	1 OF 6	REV A

TOOLS REQUIRED
1/2" WRENCH
1/2" SOCKET AND TORQUE WRENCH - OPTIONAL


ASSEMBLY COMPONENTS			
ITEMS INCLUDED	PART #	PART DESCRIPTION	QTY
TOP AND APRON ASSEMBLY	N/A	TOP ASSEMBLY (FACTORY ASSEMBLED)	1
LEGS WITH CASTERS	N/A	LEG,(OAK OR MAPLE) 2.25X2.25X29.25,HNGR BLT ATTACHED	2
LEGS FOR LEVELERS	N/A	LEG,(OAK OR MAPLE) 2.25X2.25X32.25,HNGR BLT ATTACHED	2
NUT - LEGS	100044	NUT,HEX 1/2 STEEL ZINC	4
WASHER - LEGS	100039	WASHER,SPLIT LOCK 1/2	4
LEG HARDWARE - LEVELER	520001	LEG HARDWARE KIT, 1 LEG	2
LEG BOOT	100076	LEG BOOT,BLK RUBBER	2

A	CREATED INSTRUCTIONS	09/03/25	EMR	N/A
REV.	DESCRIPTION	DATE	NAME	ECN No.
REVISION LEVEL				

ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE DECIMALS: .X = ± .030 .XX = ± .020 .XXX = ± .010 ANGLES: = ± 1° CABINET SQUARENESS = ± .0625	DWN BY:	EMR	DATE DWN:	09/03/25
	APR BY:		DATE APR:	
	APR BY:		DATE APR:	
	APR BY:		DATE APR:	
CONFIDENTIAL: ALL DESIGN RIGHTS RESERVED THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC. THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.				

			
TITLE: PERPETULAB WHEELBARROW SCIENCE TABLE TOOLS AND ASSEMBLY COMPONENTS			
DRAWING NUMBER:		SCALE:	SHEET:
P7604x34W SERIES TABLES		NTS	2 OF 6
		REVISION:	REV A

HARDWARE REQUIRED FOR STEP		
QTY	DESCRIPTION	PART NUMBER
1	TOP AND APRON ASSEMBLY	N/A

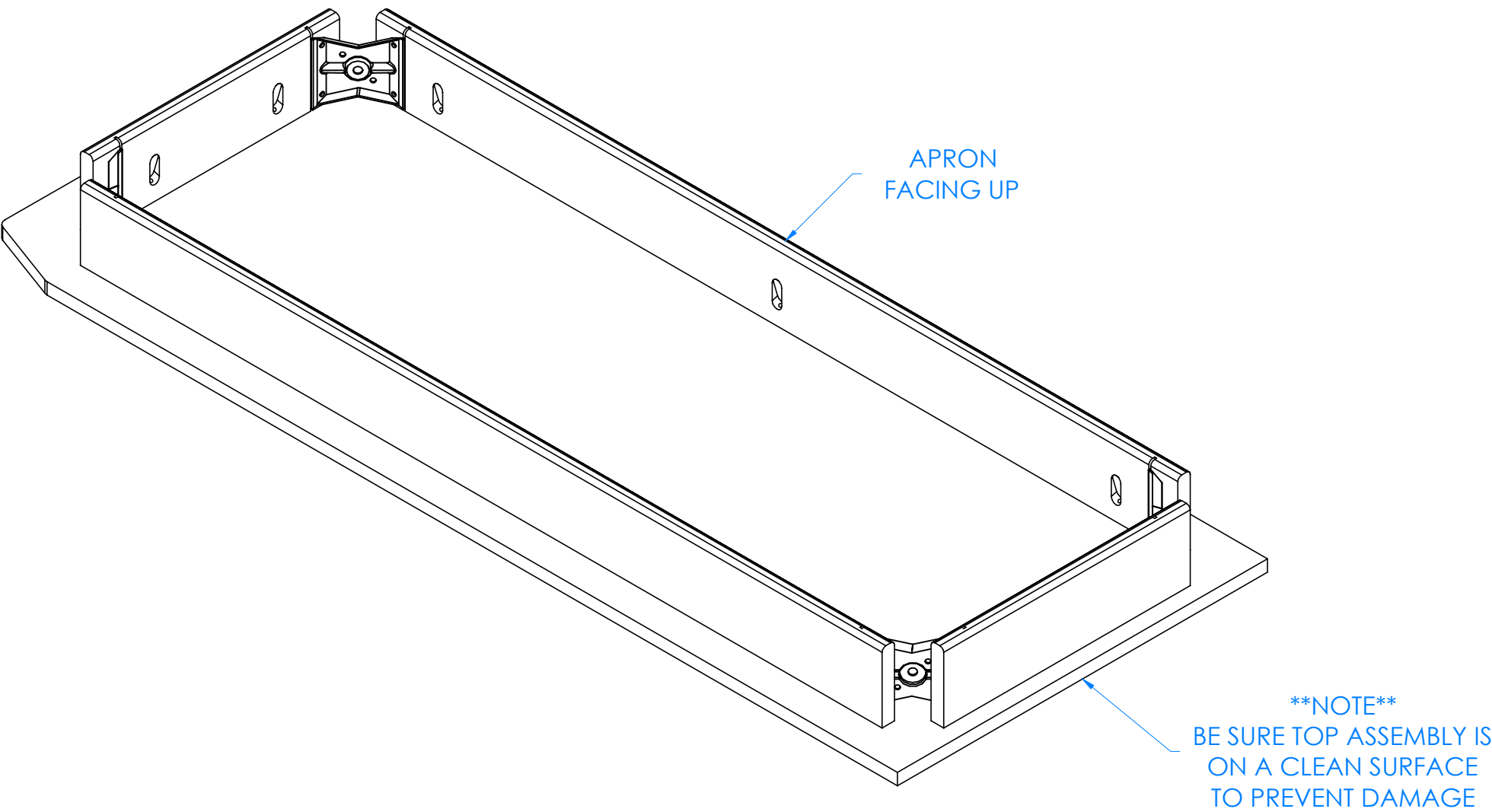

- STEP 1:
- α. LAY TOP ASSEMBLY ON FLOOR WITH APRON SIDE FACING UP.

A	CREATED INSTRUCTIONS	09/03/25	EMR	N/A
REV.	DESCRIPTION	DATE	NAME	ECN No.
REVISION LEVEL				

ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE DECIMALS: .X = ± .030 .XX = ± .020 .XXX = ± .010 ANGLES: = ± 1°0' CABINET SQUARENESS = ± .0625	DWN BY:	EMR	DATE DWN:	09/03/25
	APR BY:		DATE APR:	
	APR BY:		DATE APR:	
	APR BY:		DATE APR:	
CONFIDENTIAL: ALL DESIGN RIGHTS RESERVED THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC. THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.				



TITLE:
PERPETULAB WHEELBARROW SCIENCE TABLE
ASSEMBLY INSTRUCTIONS


DRAWING NUMBER:	SCALE:	SHEET:	REVISION:
P7604x34W SERIES TABLES	NTS	3 OF 6	REV A

STEP 2:


- a. INSERT HANGER BOLT ON ONE LEG THROUGH ONE CORNER BRACKET AND INSERT LOCK WASHER AND HEX NUT - (2A).

a.a. LEGS WITH CASTERS GO ON *OPPOSITE SIDE* OF HANDLE CUTOUT IN TOP - (2B).

a.b. LEGS WITH LEVELERS WILL GO ON SAME END WITH HANDLE CUTOUT IN TOP - (2B).
- b. SCREW ON NUT USING A 1/2" WRENCH OR 1/2" SOCKET AND TORQUE WRENCH, TIGHTEN TO A MINIMUM OF 12FT LBS BUT NOT TO EXCEED 17FT LBS.
- c. REPEAT FOR OTHER LEGS.
- d. MAKE SURE ALL LEGS ARE STRAIGHT AND ALL NUTS ARE ON TIGHT !

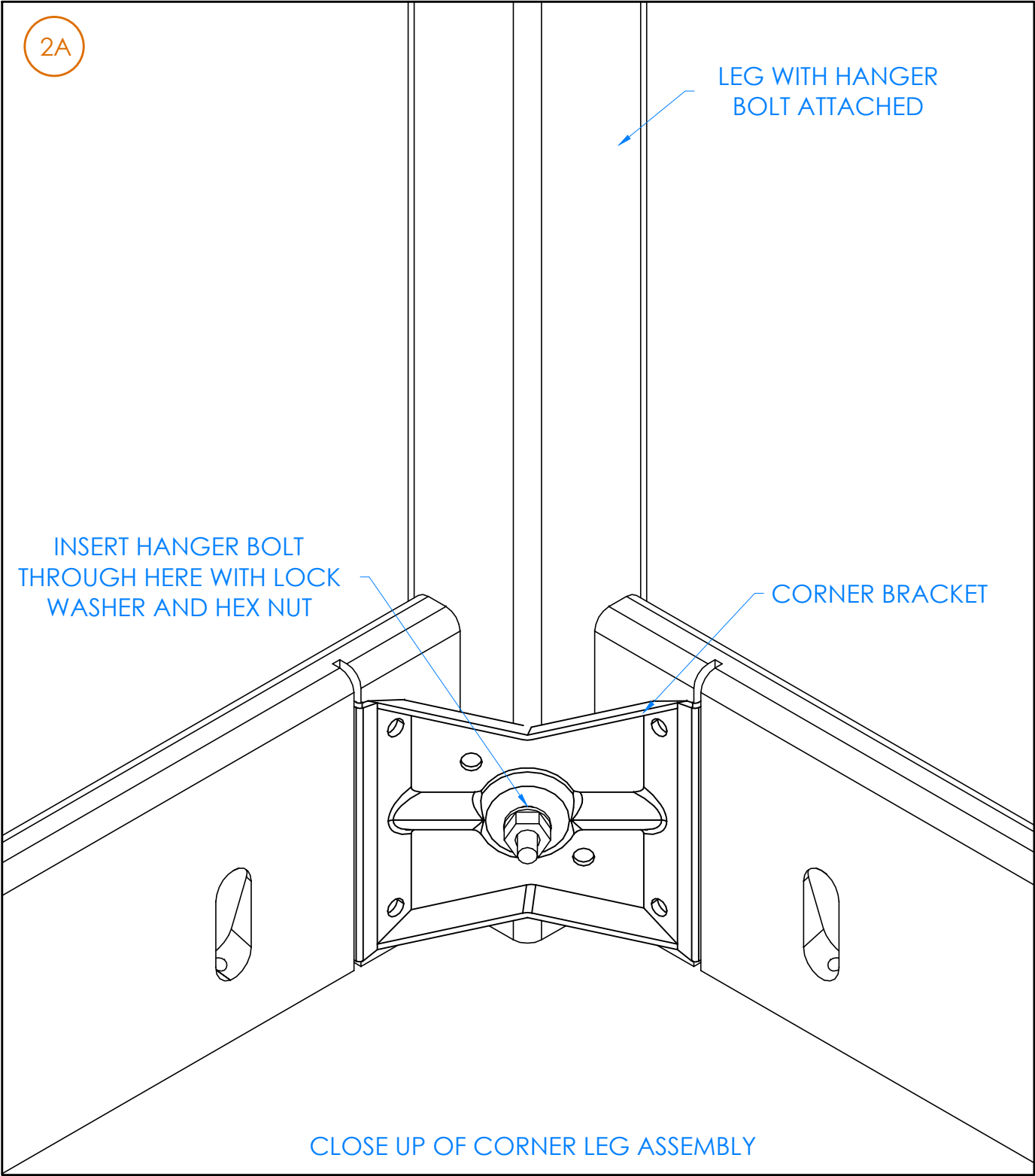

HARDWARE REQUIRED FOR STEP		
QTY	DESCRIPTION	PART NUMBER
1	1/2" WRENCH	N/A
2	LEGS WITH CASTERS	N/A
2	LEGS FOR LEVELERS	N/A
4	NUTS - LEGS	100044
4	WASHERS - LEGS	100039

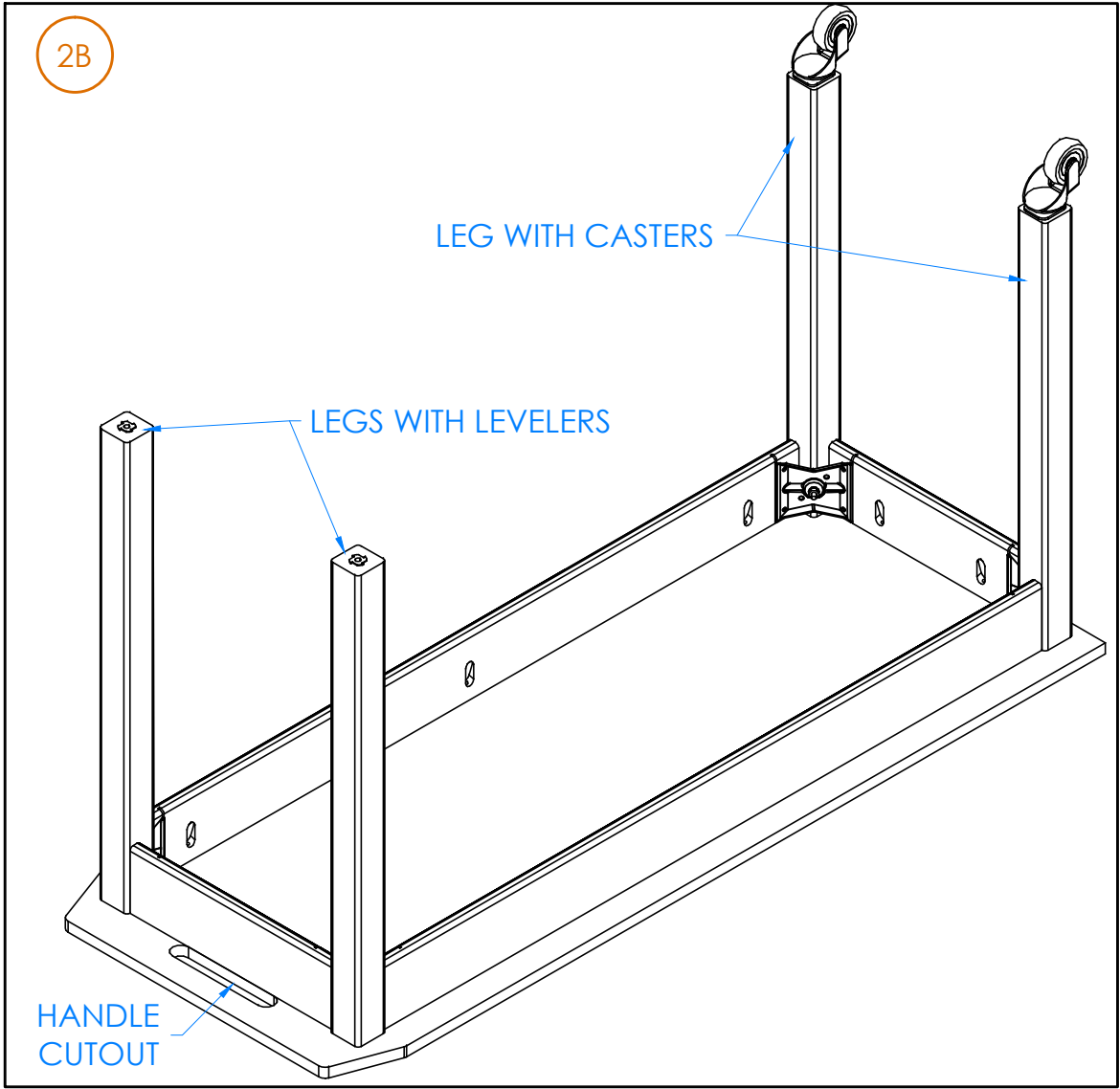

ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE
DECIMALS: .X = ± .030
.XX = ± .020
.XXX = ± .010
ANGLES: = ± 1°
CABINET SQUARENESS = ± .0625

DWN BY: EMR
APR BY:
APR BY:

DATE DWN: 09/03/25
DATE APR:
DATE APR:

CONFIDENTIAL:
ALL DESIGN RIGHTS RESERVED
THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC.

THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.



TITLE:
PERPETULAB WHEELBARROW SCIENCE TABLE
ASSEMBLY INSTRUCTIONS


DRAWING NUMBER:
P7604x34W SERIES TABLES

SCALE:
NTS

SHEET:
4 OF 6

REVISION:
REV A

A	CREATED INSTRUCTIONS	09/03/25	EMR	N/A
REV.	DESCRIPTION	DATE	NAME	ECN No.
REVISION LEVEL				

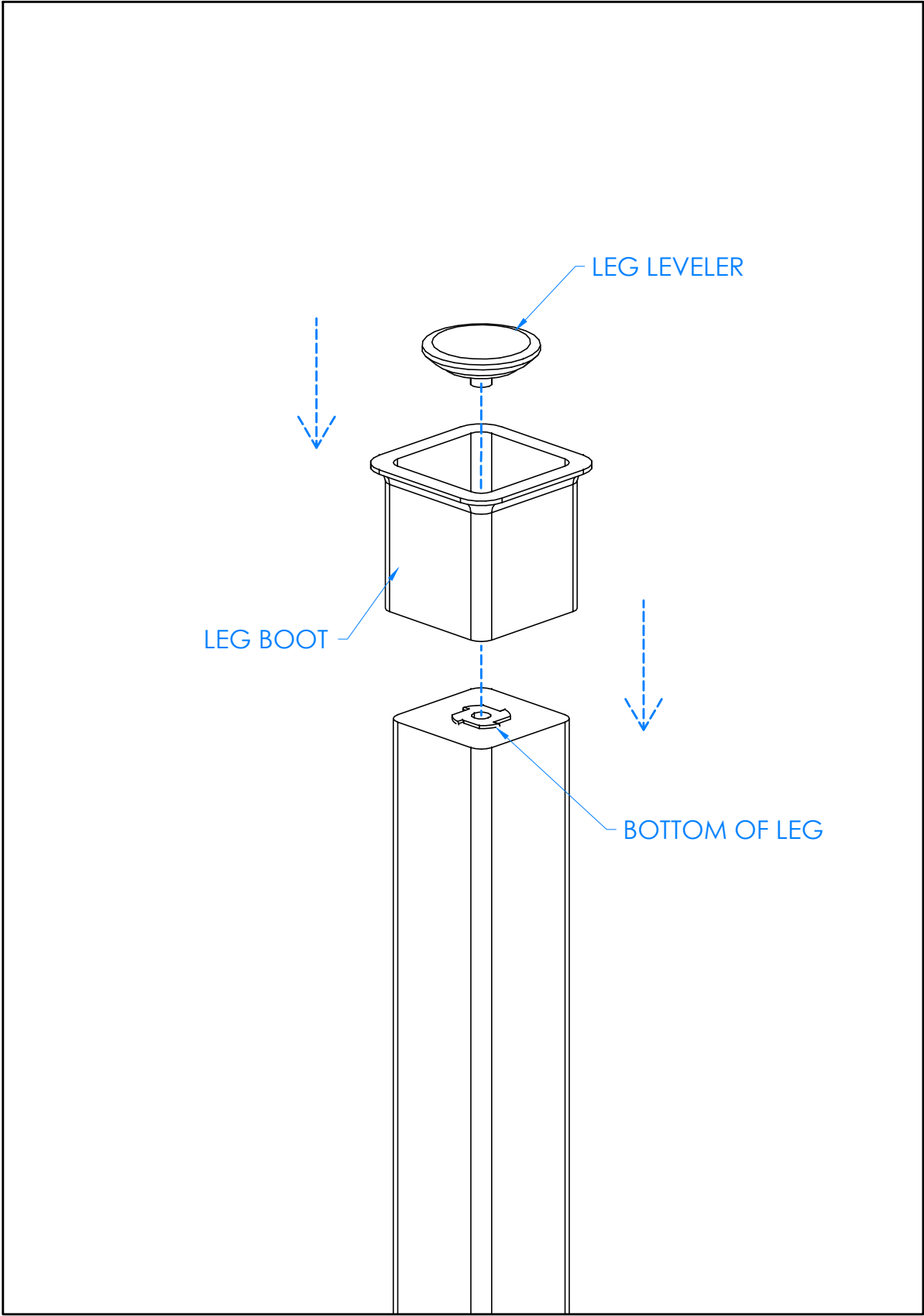

CLOSE UP OF LEG BOOT AND LEG LEVELER ASSEMBLY

HARDWARE REQUIRED FOR STEP		
QTY	DESCRIPTION	PART NUMBER
2	LEG HARDWARE - LEVELER	520001
2	LEG BOOT	100076


STEP 3:

- a. SLIDE LEG BOOTS ONTO LEGS WITH FLARED END TOWARDS BOTTOM OF LEG.
- b. THREAD LEG LEVELER INTO T-NUT LOCATED ON BOTTOM OF LEG.
- c. REPEAT FOR OTHER LEG ASSEMBLY.

ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE DECIMALS: .X = ± .030 .XX = ± .020 .XXX = ± .010 ANGLES: = ± 1°0' CABINET SQUARENESS = ± .0625	DWN BY:	EMR	DATE DWN:	09/03/25
	APR BY:		DATE APR:	
	APR BY:		DATE APR:	
	APR BY:		DATE APR:	
CONFIDENTIAL: ALL DESIGN RIGHTS RESERVED THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC.				
THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.				

			
TITLE: PERPETULAB WHEELBARROW SCIENCE TABLE ASSEMBLY INSTRUCTIONS			
DRAWING NUMBER:		SCALE:	SHEET:
P7604x34W SERIES TABLES		NTS	5 OF 6
		REVISION:	REV A

A	CREATED INSTRUCTIONS	09/03/25	EMR	N/A
REV.	DESCRIPTION	DATE	NAME	ECN No.
REVISION LEVEL				

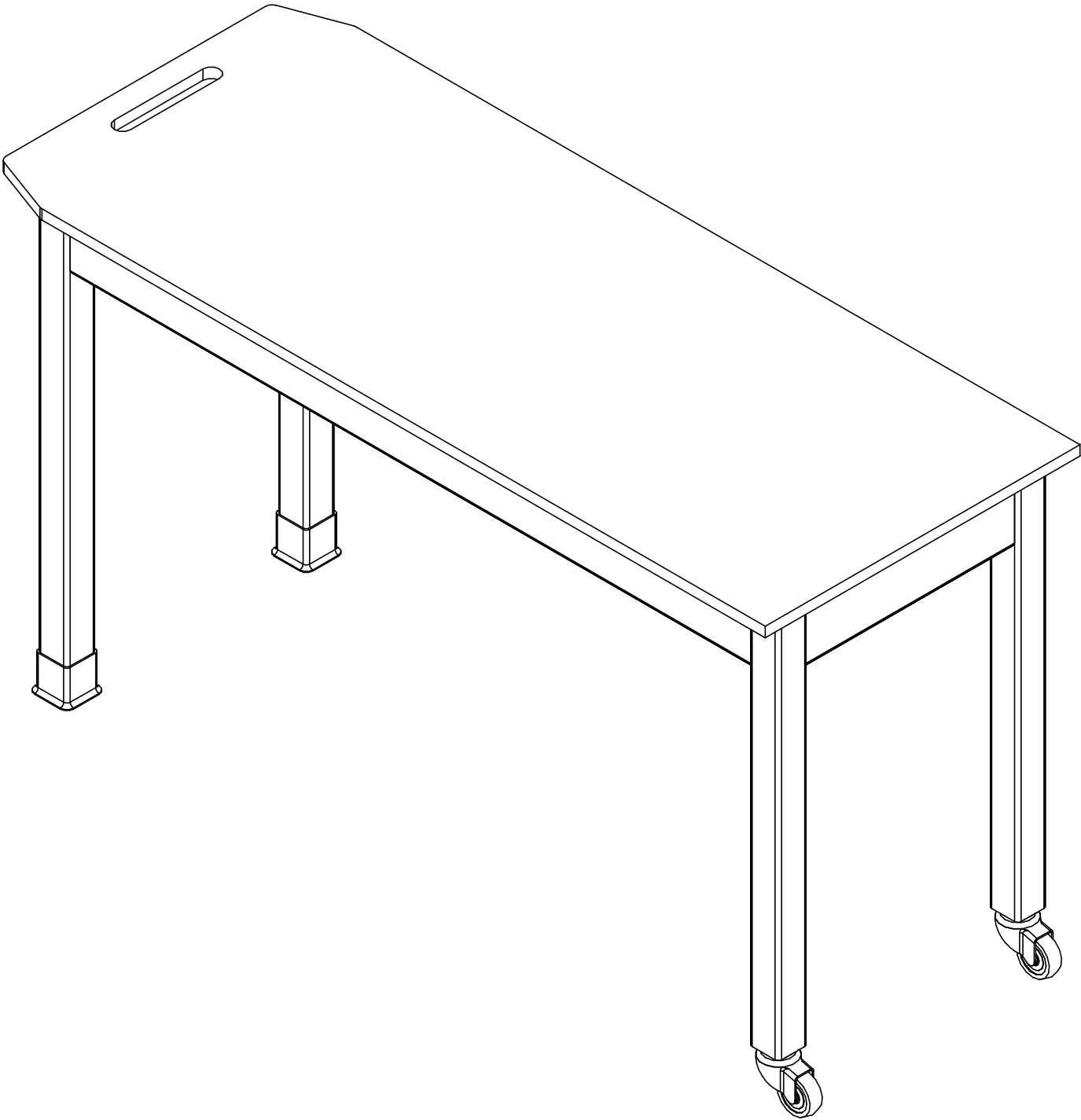

COMPLETE UNIT SHOWN

STEP 4:

- a. WITH HELP FROM MULTIPLE PEOPLE, LIFT AND TURN TABLE UPRIGHT.
- b. SET TABLE IN PLACE.
- c. ADJUST LEG LEVELERS TO LEVEL TABLE AS NEEDED.
- d. SLIDE LEG BOOTS DOWN TO FLOOR.

ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE

DECIMALS: .X = ± .030
.XX = ± .020
.XXX = ± .010

ANGLES: = ± 1°0'

CABINET SQUARENESS = ± .0625

DWN BY: EMR

APR BY:

APR BY:

DATE DWN: 09/03/25

DATE APR:

DATE APR:

CONFIDENTIAL:
ALL DESIGN RIGHTS RESERVED
THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC.

THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.



TITLE:
PERPETULAB WHEELBARROW SCIENCE TABLE
ASSEMBLY INSTRUCTIONS

DRAWING NUMBER:
P7604x34W SERIES TABLES

SCALE:
NTS

SHEET:
6 OF 6

REVISION:
REV A

A	CREATED INSTRUCTIONS	09/03/25	EMR	N/A
REV.	DESCRIPTION	DATE	NAME	ECN No.
REVISION LEVEL				